

Custom Filters without the High Costs

Learn How Semrock Optical Filters can
Help You Cut Costs, Ensure Quality,
and Optimize Your Filter Bill of Materials

Why Custom Filters DON'T have to be Costly Filters

Explore the specifications, theory, and technology behind Semrock optical filters from IDEX Health & Science and how you can reduce cost, maintain quality, and meet your filter bill-of-materials (BOM).

Optical filters are used in numerous applications, such as fluorescence microscopy, flow cytometry, point-of-care diagnostics, Raman Spectroscopy, and Optical Coherence Tomography. This blog explores the cost drivers of implementing a custom filter into an instrument to debunk the myth that custom must mean costly.

We will cover the following topics:

INTRODUCTION TO CUSTOM OVER CATALOG BENEFITS

COST DRIVERS: HIGH, MEDIUM, AND LOW

WHY PARTNER WITH SEMROCK

Written by
Toby Scrivener

Edited by
Elizabeth Bernhardt

Introduction

You've just completed your newest product design and the prototype works. The only problem? It's a massive tabletop of cluttered catalog optics with huge unused space between. By using customized optical filters designed specifically for your instrument and technique, you can:

- ▶ Reduce the size of your system
- ▶ Reduce the cost
- ▶ Reduce complexity
- ▶ Protect your supply chain

We discuss optical filter specifications, cost drivers, and reasons why you should explore custom filters as an additional cost-effective option.

Main Costs and a Collaborative Approach

At IDEX Health & Science, our mission is to assist customers through collaborative partnerships. As specification experts, we guide projects from concept to production, handle system upgrades, and accommodate non-standard packaging. Our focus is always on finding tailored solutions for your specific needs. The iterative process of optimizing specifications for application performance, while minimizing costs, involves input from various teams including customers, sales engineers, optical engineers, QA, and logistics. As shown in Figure 1, our team can divide certain specifications into high, medium, and low cost drivers. In this article, we explore these specifications and cost drivers in greater detail.

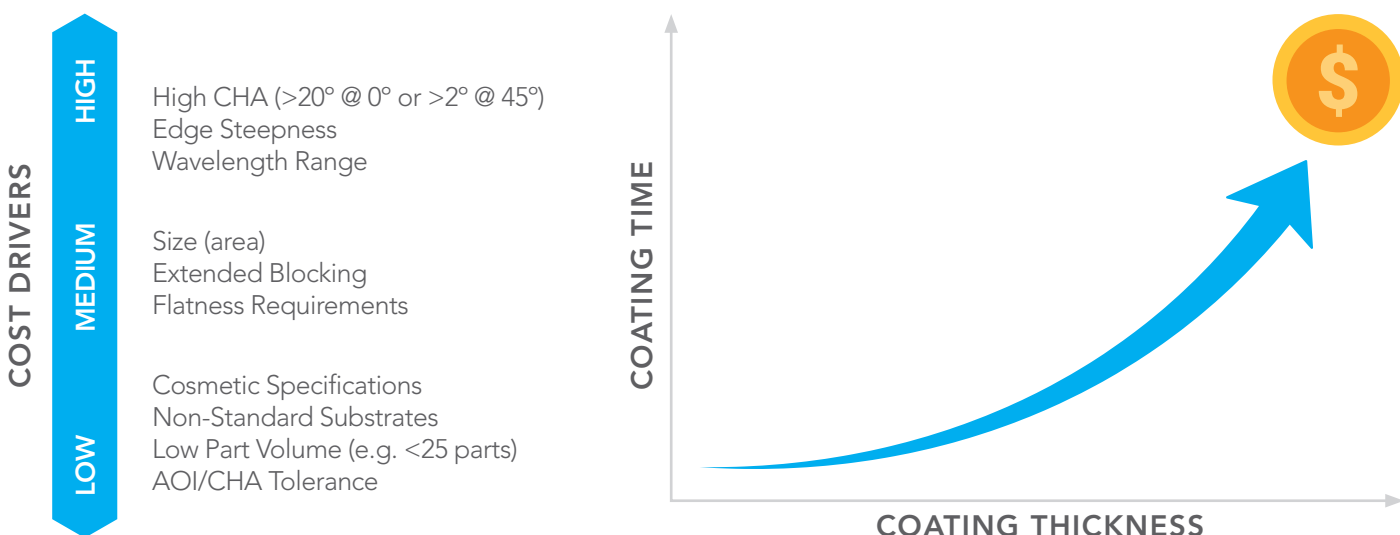


Figure 1: Optically complex responses increase coating thickness, which increases the cost.

High Cost Drivers

AOI and CHA

- › Angle of Incidence (AOI): This is the angle at which a collimated beam of light is incident on the filter, relative to the surface normal, as represented in the first two panels of Figure 2.
- › Cone Half Angle (CHA): Refers to the angular range associated with a non-collimated incident beam, as shown in the third and fourth panel of Figure 2.
- › Changing either the AOI or CHA of your incident light impacts the filter's spectral response.

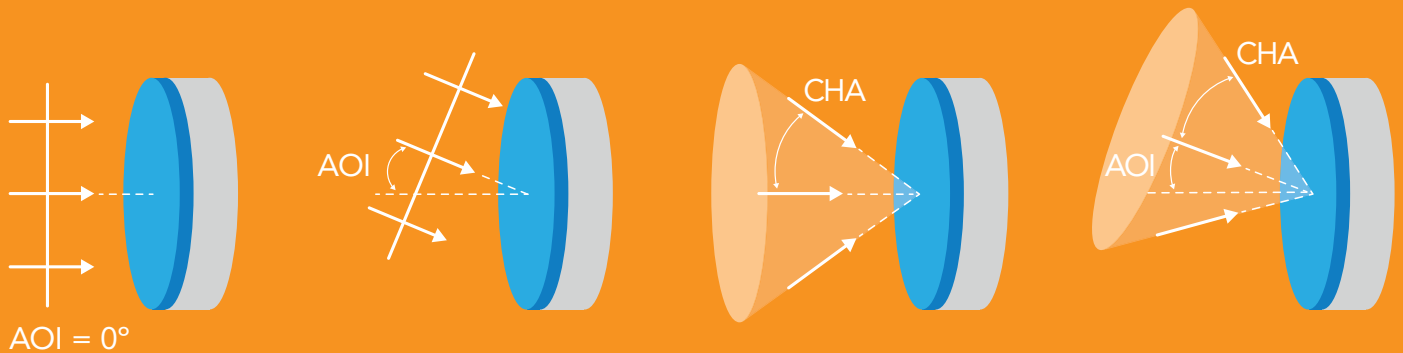


Figure 2: The difference between angle of incidence (AOI) and cone half-angle (CHA).

When discussing filter requirements, it's crucial to distinguish between Cone Half-Angle (CHA) and Angle of Incidence (AOI). Specifying an AOI range is not equivalent to specifying a filter with the same CHA value.

For instance, consider LL01-785 as a test subject. When we adjust the expected AOI of incoming light (Figure 3), the overall spectral performance of the MaxLine Laser Line filter remains relatively unchanged as the AOI increases from 0°, other than shifting to the blue as is expected for angle tuning. However, performance degrades rapidly with increasing CHA. Addressing this effect requires different design approaches.

Figure 4 illustrates the consequences of using a standard catalog filter not designed for large CHA. The desired performance (green) significantly differs from the realized performance (red). In contrast, a custom filter designed for CHA (Figure 4) shows improved out-of-band blocking and edge steepness.

Here a high CHA would be considered as anything above 20° at 0° AOI, or >2° at 45° AOI (Figure 4).

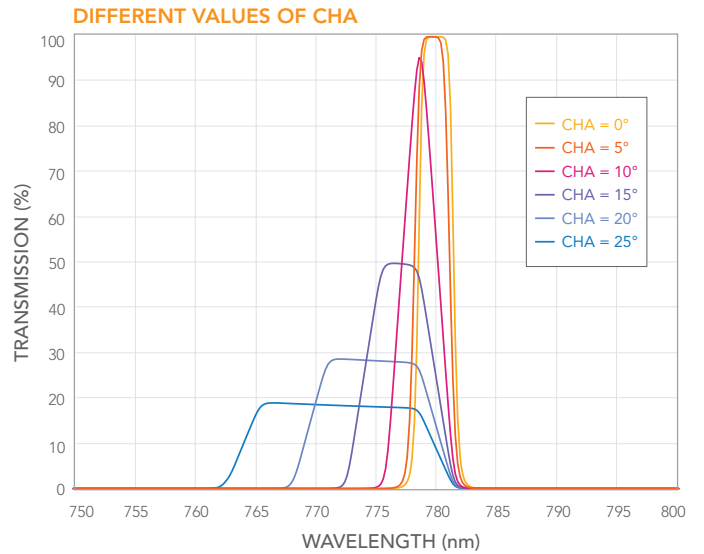
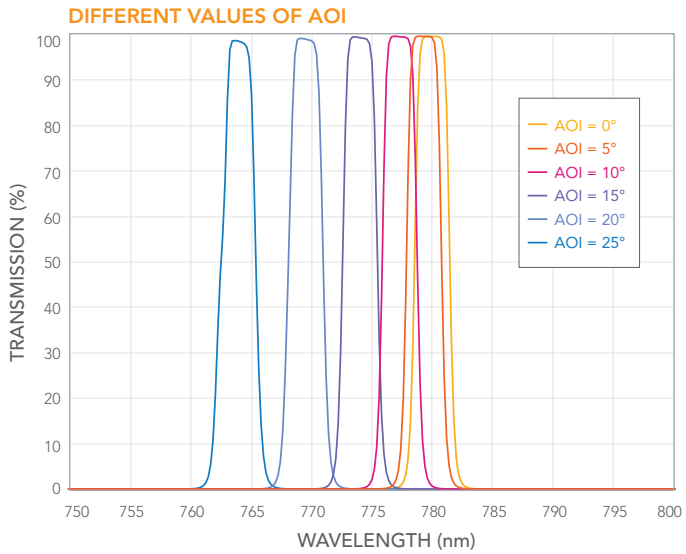


Figure 3: The effect of a varied AOI is different to a varied CHA on this MaxLine Laser Line filter.

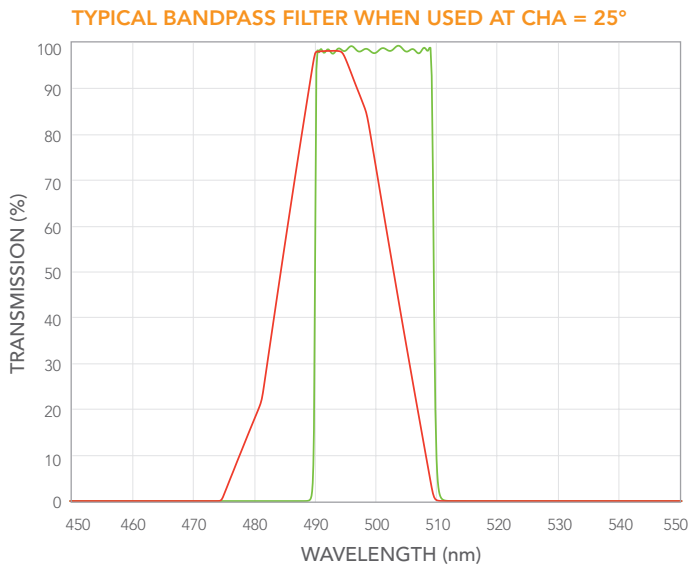


Figure 4: A standard filter without considerations to CHA shows reduced transmission performance, compared to a custom-designed filter with a high tolerance to high CHA. The green curve is CHA 0°. The red curve is CHA 25°.

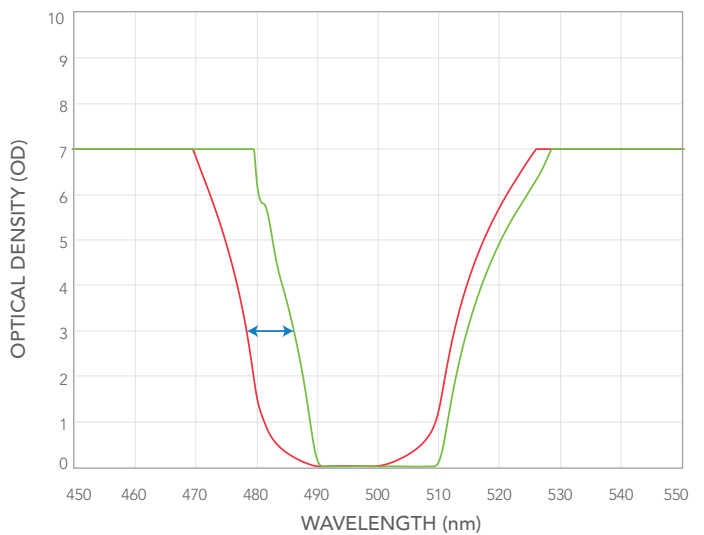
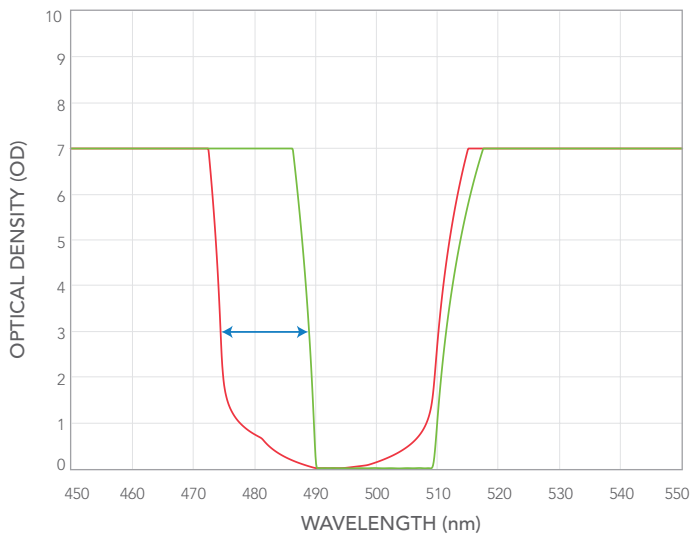


Figure 5: Out of band blocking and edge steepness is improved when CHA is compensated with a custom filter.

High Cost Drivers (Cont.)

Edge Steepness and Transition Width

Edge Steepness

Refers to the wavelength points on a filter where it transitions from blocking to transmitting.

Transition Width

This term describes the maximum distance between the laser line (where optical density, OD, is greater than 6) and the 50% transmission point.

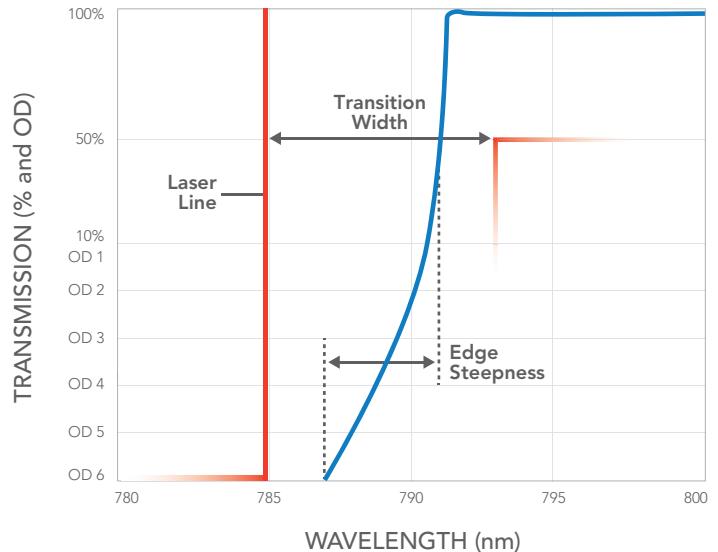


Figure 6: The difference between Transition Width and Edge Steepness.

Transition Width and Edge Steepness are related but **not interchangeable**.

Edge Steepness specifically measures the distance between where $OD > 6$ ends and the 50% transmission point.

Wavelength Range

Spectral Range

The range (bluest to reddest wavelengths, such as from 200 nm to 2000 nm) for which you need the filter designed significantly impacts costs. Moving to the extreme ends of this range can increase coating expenses.

Extreme UV

Working in the extreme UV requires unique substrates and slower deposition rates, resulting in longer production times.

IR Range

For the infrared (IR) range, we use the same materials as in the visible range, but achieving the same spectral performance requires thicker coatings. Going from 600 nm to 1200 nm doubles the required layer thickness.

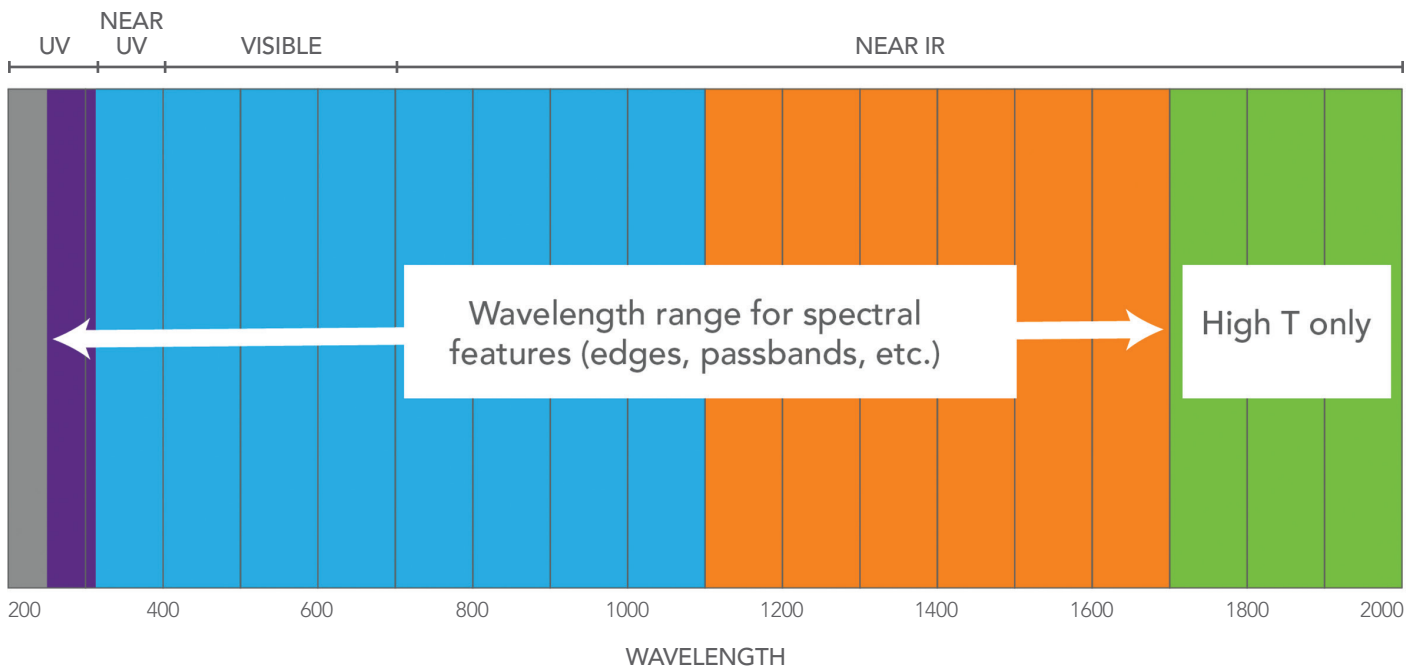


Figure 7: The spectral zones in which your features are designed have different associated costs.

Medium Cost Drivers

Flatness Requirements

Flatness refers to the deviation of a component surface from a perfectly flat reference plane. It's crucial in applications where maintaining beam shape or image quality matters. The **Reflected Wavefront Error (RWE)** measures how much a reflected wavefront deviates from a perfect one.

- At normal incidence, flatness and RWE are related by a factor of 2.
- At non-normal incidence, the relationship depends on the angle of incidence.

If you're designing a filter with strict flatness requirements, we use a stress-compensating coating. This coating balances forces to maintain flatness, but it increases cost due to longer coating chamber time. Thicker substrates can help mitigate the cost by using the number of layers needed for the stress-compensating coating.

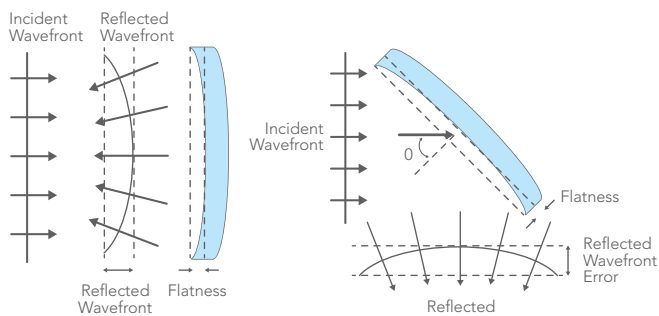


Figure 8: Examples of RWE when the optical design does not have a stress compensating optical coating.

Part Size (area)

At times, filter size can be the easiest way to reduce cost. Smaller filters generally cost less than larger ones, especially when produced in large quantities. However, very small parts (under 5 mm) pose challenges in manufacturing, handling, and packaging. Larger parts can be more complex to manufacture, and costs don't always scale linearly. During the design stage, we provide consultative advice on the optimal filter size for your application.

Extended Blocking/ Deeper Blocking

We analyze customer system requirements to recommend cost-effective blocking solutions. Often, customers ask for the highest level of blocking beyond their area of interest. By adopting a stepped approach and considering light levels and detector sensitivity, we can reduce costs significantly, as demonstrated in Figure 9.

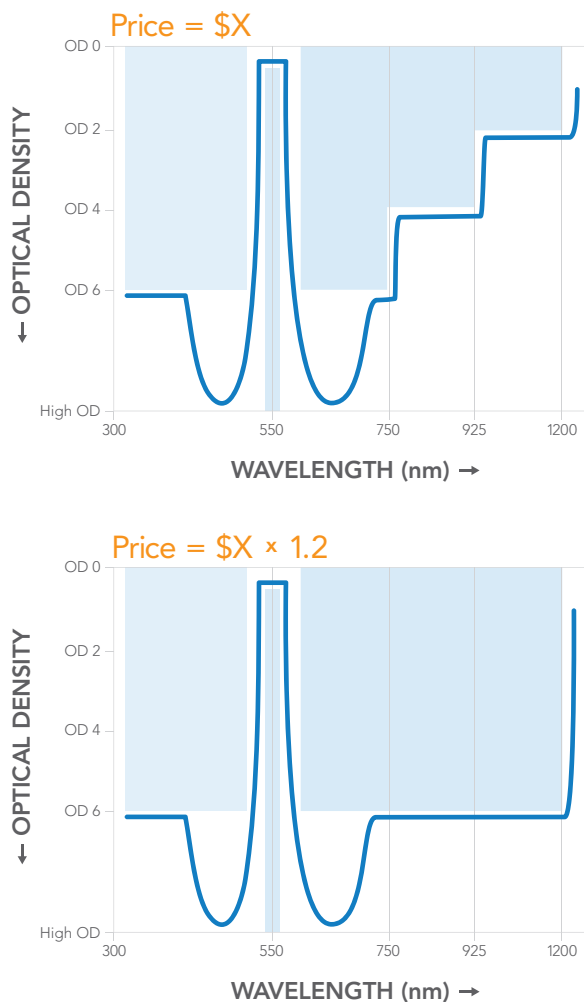


Figure 9: A balanced approach to OD (Left) has cost savings over an extended blocking approach.

Low Cost Drivers

Cosmetic Specifications

SCRATCH AND DIG

All optical coatings have cosmetic imperfections no matter how they are manufactured. There are many requirements that could be considered to be a cosmetic specification. One of the more significant specifications would be Scratch-Dig. We rely upon the definitions and inspection techniques established by the American National Standard Institute (ANSI) contained in standard ANSI/OEOSC OP1.002-2009.

Scratch and Dig (S/D):

- › Scratch: A long imperfection caused by marking or tearing on the optical surface. It has a v-groove contour and is significantly longer in dimension than it is wide. Customers specify the maximum allowable scratch width within the filter's clear aperture (CA) (e.g., a "60" designation allows a width up to 60 microns).
- › Dig: Round or irregular holes (pinholes), including grinding pits and open entrapped bubbles. Digs are characterized by their actual diameter (specified in units of 10 microns). Customers set the maximum allowable dig diameter within the CA (e.g., a "40" designation allows up to 400 micron diameter).
- › Inspection is defined by the ANSI standard. It involves using either a 40-watt incandescent lamp or two 15-watt fluorescent lamps against a black background, comparing filters by eye to a Comparison Standard at a 45° angle of incidence.
- › Only cosmetic imperfections inside the clear aperture are included in the S/D specification. More than one imperfection of each type is allowed, up to a certain number as defined by the ANSI standard.

Edge Chip and Clear Aperture

- › When parts are cut, edge chips may occur due to dislodged coating.
- › Semrock's standard specifies edge chips must lie no further than 0.5 mm from the part edge, though other specifications are possible upon request.

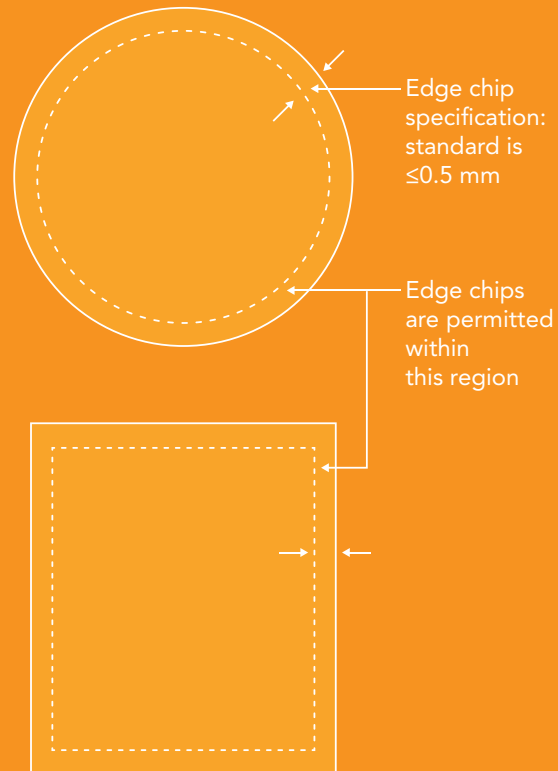


Figure 10: The CA is necessarily limited to the outer dimension minus 2x the edge chip specification.

Low Cost Drivers (Cont.)

Edge Blackening and Part Marking

Edge blackening prevents or reduces light leakage through the edges of unhoused parts. Typically done by hand for small volumes, this process can be costly. However, it may not be suitable for very small parts.

For catalog parts with standard sizes (25 mm \varnothing or 12.5 mm \varnothing), an aluminum ring housing can be used. These rings can be laser marked with

custom text and an arrow indicating the direction of light propagation. Alternatively, for parts that don't fit into a housing ring, handwritten text or a caret can serve the same purpose.

Tighter cosmetic specifications impact cost by reducing the yield but are available upon request.

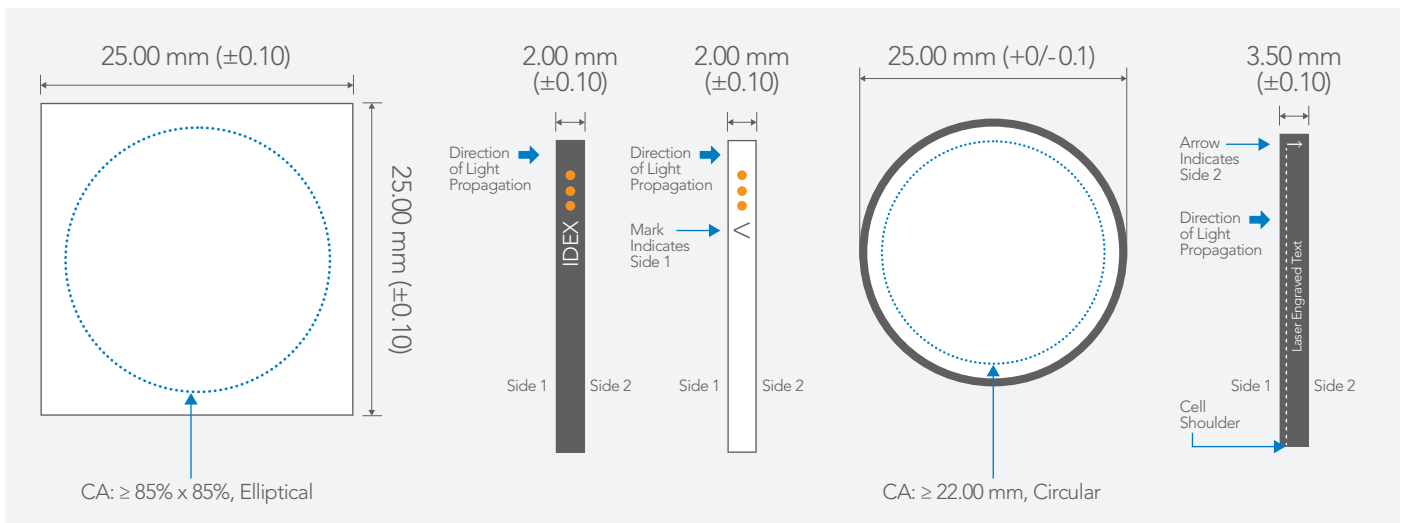



Figure 11. The options for part marking are numerous.

Non-Standard Substrates

If your optical filter needs a non-standard substrate, additional substrate polishing may be required to balance the polishing table. Without this, meeting target specifications is unlikely, leading to increased costs and lead time.



Single Substrate Can Lead to Unbalanced Table
Substrate won't meet specifications.

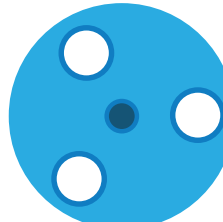

Typically 3 to 5 Substrates Are Polished
Balanced table results in specifications being met.

Figure 12. Multiple substrates are required to balance the polishing process.



AR (2nd Side Coating)

Applying an AR coating to one side of a substrate adds to the cost of creating an optical filter. However, since these coatings are thinner than the primary optical coating, they take less time to apply, making them a low-cost driver.

Other Considerations

We can consider other important qualities and specifications, such as GDD Compensation, Zero Pixel Shift, and Polarization specifications. Discuss any significant considerations for your application with your sales representative.

Why Partner with Semrock?

Semrock Optical Filters from IDEX Health & Science set industry standards by combining advanced ion-beam-sputtering systems with proprietary deposition control technology, predictive algorithms, and volume manufacturing. This results in unmatched performance for biotech and analytical instrumentation.

Customer Satisfaction

Our team continuously improves supply chain management and manufacturing processes, ensuring rigorous documentation, revision control, and traceability. We maintain inventory levels to meet delivery goals.

Flexible Delivery

We offer filters to meet demanding specifications, available through blanket orders, purchasing contracts, or just-in-time delivery. Our products cater to a wide range of applications, including:

- › Fluorescence filters
- › Raman spectroscopy filters
- › Tunable bandpass filters
- › Deep notch filters
- › Laser-line filters
- › Laser diode clean-up filters
- › Filters for combining or separating laser beams
- › Filters for isolating mercury lamp lines
- › Polarization filters
- › Dispersion controlled coating designs
- › Laser mirrors

Custom Solutions

Our core strengths include:

- › Wavelength functionality from 230 nm to 2000 nm
- › Production capacity of tens of thousands of parts per month
- › Proprietary spectral measurement systems such as KolaDeep to measure to OD8+.
- › Sizes from 1 mm² on 0.5 mm thick glass to 200 mm
- › Spectrally complex custom designs (e.g., LED optimized filters)
- › Custom sizing (round, square, rectangular)
- › On-filter laser engraving for easy identification and storage



Custom Filters

Why choose custom optics with Semrock?

Semrock catalog parts are world-renowned for their performance, and we offer custom-sized parts as a standard service. So, why switch to custom parts? Here are four reasons:

Reduce System Size

Custom optics designed to your specifications can fit more compactly, allowing you to create instruments with a smaller footprint or fit more optics into the same space.

Reduce Costs

By consulting on required specifications and considering cost drivers, we can make custom filters as affordable as standard catalog parts.

Reduce Complexity

Custom filters tailored to your optical requirements can reduce the number of filters needed, improve performance, and create an optimized beam path.

Protect Your Supply Chain

Custom filters with long-term contracts ensure your parts always meet specifications, avoiding issues from changes to catalog parts. We consider your needs when making any changes in our coating facility.

Whether from the first lot or the last, our filters maintain consistent spectral properties. OEM customers can rely on a secure and dependable supply line, as shown by the reproducible performance of the same filter from 40 different coating runs in Figure 13.

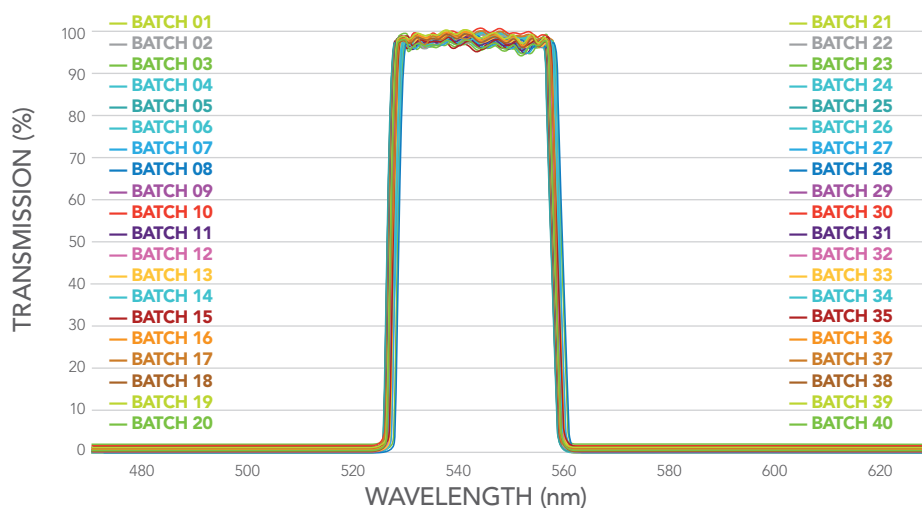


Figure 13. Forty coating runs in different coating chambers and facility locations demonstrating reproducibility and reliability.




If you're looking to:

- › REDUCE THE SIZE OF YOUR SYSTEM
- › CUT COSTS WITHOUT COMPROMISING QUALITY
- › SIMPLIFY YOUR OPTICAL DESIGN
- › SECURE YOUR SUPPLY CHAIN WITH RELIABLE, CUSTOM SOLUTIONS

Semrock is here to help you achieve your goals. With our expert team and advanced technology, we'll collaborate with you to design filters that meet your exact specifications.

Ready to explore custom or standard filter options?

Contact us today at
www.idex-hs.com/custom-optics

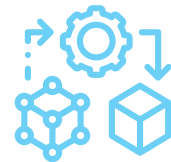



For ordering, technical support,
and contact information please visit
www.idex-hs.com/semrock